

Алматы (727)345-47-04
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Россия +7(495)268-04-70

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Казахстан +7(727)345-47-04

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Беларусь +(375)257-127-884

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саранск (8342)22-96-24
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Узбекистан +998(71)205-18-59

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Киргизия +996(312)96-26-47

<https://gonotec.nt-rt.ru> || gcd@nt-rt.ru

Осмометры для определения точки замерзания серии OSMOMAT 3000

OSMOMAT® 3000 Series

Freezing Point Osmometer



AREAS OF APPLICATION

Medical

- ▶ Clinical and diagnostic laboratories
- ▶ Neonatology and pediatrics
- ▶ Intensive care
- ▶ Urology and nephrology
- ▶ Transfusion medicine and dialysis
- ▶ Gynecology and IVF
- ▶ Medical research


Pharmacy and biotechnology

- ▶ Quality control of pharmaceutical formulations
- ▶ Bioprocess engineering
- ▶ Drug research
- ▶ Synthetic biology

Other areas of application

- ▶ Veterinary medicine
- ▶ Various research areas
- ▶ Food and beverage industry and analytics
- ▶ Agriculture
- ▶ Consumer goods industry
- ▶ Environmental analytics
- ▶ Chemical industry
- ▶ Forensic medicine

The perfect laboratory companions to determine the osmolality of aqueous solutions

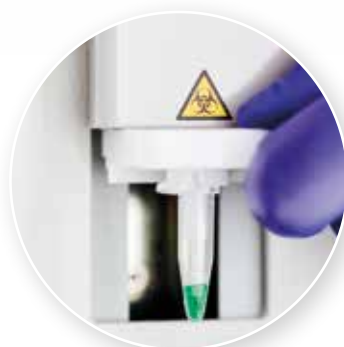

Software for data transfer

For serial data transfer to a PC or information system (like LIS or HIS) a corresponding receiving software is needed on the target system. GONOTEC does not supply such software solutions and does not support software products from other manufacturers.

APPLICATION

Freezing point osmometers are used in numerous areas of application, such as medical, pharmaceutical, biotechnology, food and beverage, and chemical industries.

In the human body osmotic processes and osmoregulation play an important role: a disturbance of the osmotic equilibrium can lead to numerous health impairments. In many cases, osmometers are used in the medical and pharmaceutical industries to determine the osmolality of blood or urine samples or pharmaceutical formulations by freezing point osmometry.

The **Gonotec® OSMOMAT® 3000 series** have been specially designed for routine **measurements in the medical and pharmaceutical fields** but due to their robustness, precision and ease of use the devices are the perfect choice for many others areas as well.

The **OSMOMAT® 3000** determines the total osmolality of aqueous solutions, requires extremely small sample amounts, and allows serial measurements in the shortest possible time.

MEASUREMENT TECHNOLOGY

Osmometry is an analytical measuring method for determining the osmotic value / pressure (often referred to as simply the osmolality) of a sample.

The osmolality is defined as the concentration of all dissolved – and thus osmotically effective – particles in a solution based on 1 kilogram of solvent. **The unit of osmolality is Osm/kg or Osmol/kg.**

The freezing point of a sample changes depending on the concentration of dissolved substances.

Pure water has a freezing point of 0 °C. The solution of one or several substances in water leads to lowering of the freezing point. A solution with an osmotic value of 1 Osmol/kg has a freezing point of -1.858 °C.

Through the linear correlation between the freezing point of a sample and its osmolality, **freezing point osmometry provides high-precision analytics.**

SPECIFICATIONS OSMOMAT® 3000 Series

Models	3000 <i>basic</i> , 3000, 3000 M, 3000 D, 3000 D-M
Display	5.7" LCD touch screen
Weight	6.5 kg (14.3 lbs.)
Dimensions (WxHxD)	205 mm x 360 mm x 220 mm (8.1" x 14.2" x 8.7")
Cooling	2 separate peltier elements / heat dissipation through active ventilation
Sample volume	50 µL / 15 µL for models 3000 M und D-M
Measurement	Measurement of single samples or batch processing (only single sample measurement for model 3000 <i>basic</i>)
Measurement time	~ 60 seconds
Resolution	1 mOsmol/kg H ₂ O
Units	mOsmol/kg, Osmol/kg, °C
Measurement range	0 bis 3000 mOsmol/kg H ₂ O
Reproducibility Models 3000 <i>basic</i> , 3000 und 3000 D	≤ 2 mOsmol/kg (SD) [0 bis 400] mOsmol/kg ≤ 0,5 % (CV) [400 bis 1500] mOsmol/kg ≤ 1 % (CV) [1500 bis 3000] mOsmol/kg
Reproducibility Models 3000 M und D-M	≤ 4 mOsmol/kg (SD) [0 bis 400] mOsmol/kg ≤ 1 % (CV) [400 bis 1500] mOsmol/kg ≤ 2 % (CV) [1500 bis 3000] mOsmol/kg
Calibration	2 point calibration, 3 point calibration
Linearity	Deviation less than ±1 % in the calibrated range
Ambient temperature	10 °C to 35 °C
Power supply	100 – 240VAC, 50 / 60 Hz, 80 VA
Interfaces	RS-232, USB (not for model 3000 <i>basic</i>)
Output formats	CSV, XML (not for model 3000 <i>basic</i>)
Printer (optional)	Graphical dot matrix printer for date, time and sample information for each measurement (for models 3000 D and 3000 D-M)
Printer paper	Plain paper, 43 mm (1.7") wide
Ribbon	Endless ink ribbon cartridge, replaceable
Error messages	Printed in plain text
Software laguages	German, English, Spanish, French, Portuguese, Chinese

Legend

SD: Standard deviation, CV: coefficient of variation (= relative SD)

	Item number
OSMOMAT® 3000 <i>basic</i>	# 32.B
OSMOMAT® 3000	# 32.00000
OSMOMAT® 3000 M	# 32.02000
OSMOMAT® 3000 D	# 32.10000
OSMOMAT® 3000 D-M	# 32.12000

OPTIONAL ACCESSORIES

	Item number
Handheld barcode scanner with power supply, connection cable and manual	35.9.2000

ACCESSORIES AND SUPPLIES

Accessories	Item number
250V Power Cord - Europe Plug Typ E+F (CEE 7/7)	20.9.0100
RS-232 data cable	20.9.0165
USB cable	20.9.0166
Adjustment tool	30.2.0030
Ampoule opener	30.9.1050
Bellow (pasteur pipette)	30.9.0030

Supplies	Item number
Calibration Standard 100 mOsmol/kg NaCl/H ₂ O, 10 x 1 ml	30.9.0100
Calibration Standard 300 mOsmol/kg NaCl/H ₂ O, 10 x 1 ml	30.9.0020
Calibration Standard 500 mOsmol/kg NaCl/H ₂ O, 10 x 1 ml	30.9.0500
Calibration Standard 850 mOsmol/kg NaCl/H ₂ O, 10 x 1 ml	30.9.0850
Calibration Standard 2000 mOsmol/kg NaCl/H ₂ O, 10 x 1 ml	30.9.2000
Reference Solution OSMOREF® 290 mOsmol/kg NaCl/H ₂ O, 10 x 1 ml	30.9.0290
Printer paper, OSMOMAT® 3000 D und D-M / 8 rolls	30.9.1010
Endless ink ribbon cartridge, OSMOMAT® 3000 D and D-M	30.9.1020
Measuring vessels / 1,000 pcs.	30.9.0010

STANDARD ACCESSORIES, *included*



- 1 Power cord
- 2 RS-232 data cable (not for model 3000*basic*)
- 3 USB cable for connection to PC or Laptop (not for model 3000*basic*)
- 4 Measuring vessels, 100 pcs.
- 5 2 spare fuses T 1.6A (HBC 1500A)
- 6 Adjustment tool
- 7 Calibration Standard 300 mOsmol/kg, 10 x 1 ml
- 8 Calibration Standard 850 mOsmol/kg, 10 x 1 ml
- 9 Ampoule opener
- 10 Bellow (pasteur pipette)
- 11 User guide (on compact disc (CD))

MEASUREMENT METHOD: Freezing Point Osmometer



Position measuring vessel on the thermistor probe.


Sample is moved into the lower cooling system and cooled to less than 0 °C by a microprocessor-controlled peltier element.

Injection of an ice crystal by cooled triggering needle (cryst-needle).

Crystallization of the sample material.



At the beginning of the measurement, the sample is supercooled in the lower cooling system. The freezing process of the sample material is released in a controlled manner by injecting an ice crystal using the release/triggering needle. The crystallization process leads to the release of thermal energy and the temperature of the sample increases until a plateau phase or a maximum is reached, which depicts the actual freezing point of the sample. At this point, the temperature is measured with an accuracy of 0.001 °C using a high-precision temperature sensor.

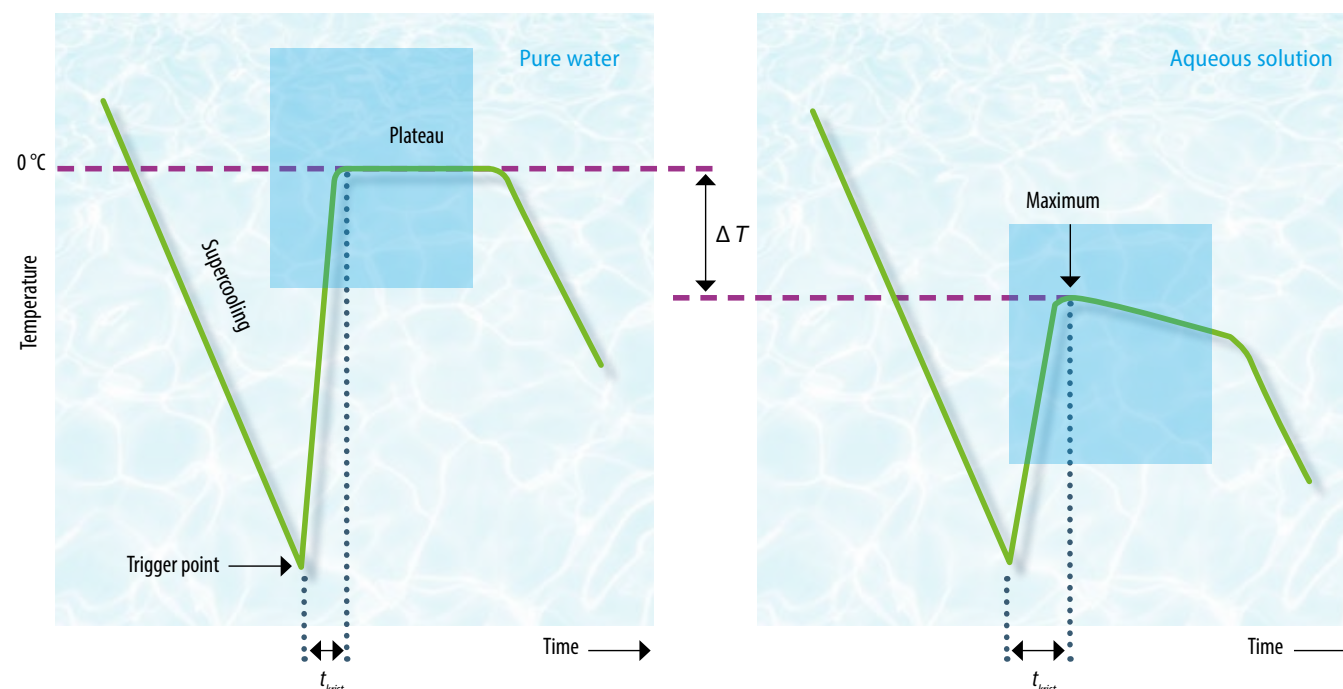

OSMOMAT® 3000basic

The Minimalist Model

- ▶ Sample volume: 50 µL
- ▶ Easy handling and maintenance
- ▶ Comfortable calibration routine
- ▶ No connectivity
- ▶ Delivery including standard accessories



OSMOMAT® 3000 / 3000 M

The Standard Models

- ▶ Sample volume: 50 µL (optional 15 µL / for model 3000 M)
- ▶ Easy handling and maintenance
- ▶ Comfortable calibration routine
- ▶ Connectivity for data transfer to PC or laptop
- ▶ Optional barcode reader
- ▶ Delivery including standard accessories and qualification document (IQ/OQ/PQ)



OSMOMAT® 3000 D / 3000 D-M

The Premium Models with Printer

- ▶ Sample volume: 50 µL (optional 15 µL / for model 3000 D-M)
- ▶ Integrated printer (delivery includes 1 ink ribbon cartridge and 1 printer paper roll)
- ▶ Easy handling and maintenance / comfortable calibration routine
- ▶ Connectivity for data transfer to PC or laptop
- ▶ Optional barcode reader
- ▶ Delivery including standard accessories and qualification document (IQ/OQ/PQ)



Your advantages at GONOTEC – AN ELITECHGROUP COMPANY

OSMOMAT® 3000 Product Series

- ▶ Robust, precise, reliable, fast
- ▶ Small sample volume
- ▶ Easy control via the integrated touch screen display
- ▶ Step-by-step guidance through all measuring functions and setting options
- ▶ 2 or 3 point calibration
- ▶ Automatic and safe calibration with the Gonotec Calibration Standards
- ▶ Data transfer to PC or laptop via RS-232 or USB (not for model 3000basic)
- ▶ Quality made in Germany
- ▶ Over 40 years of experience in precision measurement technology and osmometry
- ▶ Comprehensive advice and reliable service from our experts

Алматы (727)345-47-04
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саранск (8342)22-96-24
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(727)345-47-04

Беларусь +(375)257-127-884

Узбекистан +998(71)205-18-59

Киргизия +996(312)96-26-47

<https://gonotec.nt-rt.ru> || gcd@nt-rt.ru